# Data sheet

## **Firefly Luciferase Assay Kit**


### Cat. No: CA130 (100 assays) Cat. No: CA135 (1000 assays)

### Introduction

Luciferase from the firefly (Photinus pyralis) is the most commonly used photoprotein in molecular biology studies and has been used as a reporter for studying gene regulation and function in transformed cell lines in culture.

Firefly luciferase has an apparent molecular weight of 62 kDa, which is active as a monomer and does not require subsequent processing for its activity. The enzyme catalyzes the oxidation of reduced luciferin in the presence of ATP-Mg2+ and oxygen to generate CO<sub>2</sub>, AMP, PPi, oxyluciferin, and produces a flash of light that is proportional to the quantity of luciferase in the reaction mixture.

The Luciferase Assay Substrate includes coenzyme A, ATP, and luciferin. Including coenzyme A in the reaction enhances the sensitivity of the assay and provides a sustained light reaction (half-life >5 minutes). This eliminates the need for automated luminometer injection of substrate and allows analysis by photographic film or scintillation counting.

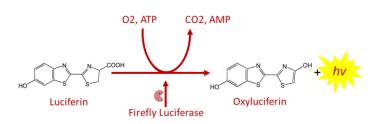

The Canvax Firefly Luciferase Assay Kit is designed for simple and efficient quantitation of firefly luciferase reporter enzyme activity from cultured cells with high sensitivity and linearity.

### **Kit Contents**

	CA130	CA135
Luciferase Assay Substrate	10 ml	100 ml
5X Cell Lysis Buffer*	4ml	20 ml
Luciferase (control)	10 µg	50 µg

\* Prepare 1X Cell Lysis Buffer by adding 4 volumes of water to 1 volume of 5X Cell Lysis Buffer.

### Storage condition

The kit is shipped on gel pack. Store all kit components at -20°C or below, protected from light. The kit is stable at -20°C for three months and at -70°C for up to 1 year from date of receipt.

Avoid repeated freeze-thaw cycles.

### **Advantages/Features:**

- Easy to use.
- Rapid: results within 15 minutes.
- Sensitivity
- Accuracy
- Ideal for high throughput assays.

### (continued on reverse side)

Distributed by:



Tallaght Business Park Whitestown, Dublin 24, Ireland D24 RFK3

Tel: (01) 4523432 Fax: (01) 4523967 Web: www.labunlimited.com Quatro House, Frimley Road, Camberley, United Kingdom **GU16 7ER** 

Tel: 08452 30 40 30 Fax: 08452 30 50 30 E-mail: info@labunlimited.com E-mail: info@labunlimited.co.uk Web: www.labunlimited.co.uk



### **Assay Procedure**

Note: The Luciferase Substrate Solution and samples should be at ambient temperature prior to performing a luciferase assav.

The reagent has been validated in a **96-well** format. Other format will require scaling and optimization by the end-user.

### A. Preparation of Cell Extract

canvax

- Aspirate the growth medium from cells. Wash cells with PBS. 1.
- 2. Add 50 µl of Lysis Buffer 1X per well of a **96-well plate**. Incubate at room temperature for 10-15 minutes.

### **B. Standard curve**

3. Produce a standard curve of light emission versus enzyme concentration.

To produce a standard curve of light units versus relative enzyme concentration, make serial dilutions of recombinant luciferase (standard) in 1X lysis buffer supplemented with 1mg/ml BSA. The addition of BSA is necessary to ensure that luciferase is not lost from solution by adsorption.

a. Dilute Luciferase to 1000 ng/ml in lysis buffer.

- b. Make serial dilutions to get concentration of 5, 2.5, 1.25, 0.625, 0.312, 0.156 ng/ml and a blank control.
- c. Add 20 µl of serial dilutions to the wells of a 96-Well Solid White Plate. (For a 384- Well Solid White Plate add 10  $\mu$ l of each concentration to the Well)

### **C. Assay Protocol**

- Transfer 20 µl of cell extracts containing Luciferase to the corresponding well of a 96-Well Solid Plate (white) with 4. lid.
- Add 100 µl/well of substrate solution in each well of cell extract (or luciferase standard). Shake 30 seconds the plate 5. to get homogenize the reaction.

(For **384-well plates**, add 25  $\mu$ l of the Luciferase substrate solution to 10  $\mu$ l of the cell lysate or luciferase standard)

- 6. Program the luminometer for the appropriate delay and measurement times.
- 7. Measure the light produced.

### D. Plot the Standard Curve

### E. Determine the Sample Concentration

- 8. Use the blank control to eliminate the background. Measure the samples.
- Extrapolate the luminescence data (light units) to Luciferase standard curve to calculate de concentration of the 9. enzyme in the cell extracts.

### **PRODUCT USE LIMITATION**

This product is developed, designed and sold exclusively for research use only. Not for diagnostic or therapeutic use. Please refer to www.canvaxbiotech.com for the Material Safety Data Sheet of the product.

### Distributed by:



Tallaght Business Park Whitestown, Dublin 24, Ireland D24 RFK3

Tel: (01) 4523432 Fax: (01) 4523967 Web: www.labunlimited.com

Quatro House, Frimley Road, Camberley, United Kingdom GU16 7ER

E-mail: info@labunlimited.com E-mail: info@labunlimited.co.uk

Tel: 08452 30 40 30 Fax: 08452 30 50 30 Web: www.labunlimited.co.u

