


## Super Clean Products

**Smart Gas Purification Products – Super Clean Gas Purifiers** 

ľ

# SGT technology creates Simply Smart Products

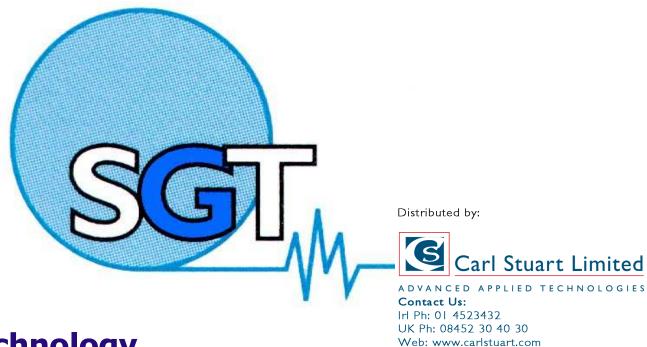

# **Scientific Glass Technology**

Middelburg, The Netherlands - Baden-Baden, Germany - Singapore

03.2007

Email: info@carlstuart.com



## Super Clean Products

### **Smart Gas Purification Products – Super Clean Gas Purifiers**

2



#### **Rear end connectors:**

- Brass or SS
- 1/4" or 1/8"

#### **INSTRUCTIONS** how to replace a CONNECTOR:

- Remove the connector, which should be replaced, from the base plate
- 2. Remove copper O-ring from its seat in the hole
- 3. Insert a new copper O-ring in the seat in the hole
- 4. Assemble the new connector into the base plate with a torque of 30 Nm
- Perform a leak test on the new connector and the base plate with a leak detector, Snoop etc.

#### Remark:

Since SGT is using a <u>plan parallel thread</u> for Connectors and Base Plates, it is POSSIBLE to replace connectors; some other manufacturers are using PNT thread, where it is IMPOSSIBLE to replace the connector!!