Data sheet

Carbenicillin disodium salt

Cat. No: AB007 Cat. No: AB008

Introduction

Carbenicillin Disodium Salt is a semi-synthetic penicillin antibiotic with a relatively greater spectrum of activity against Gram-negative bacteria. As penicillin, is susceptible to β-lactamase, which hydrolyzes the β-lactam ring. Ampicillin is widely used in cell culture as a selective Inhibits bacterial cell-wall synthesis inactivating transpeptidases on the inner surface of the bacterial cell membrane.

α-Carboxybenzylpenicillin disodium salt

Technical specification

✓ CAS Number: 4800-94-6

MDL: MFCD00077683

✓ Formula: C₁₇H₁₆N₂Na₂O₆S

Formula Weight: 422.36 g/mol

✓ Soluble in water (50 mg/ml)

Product Size

| | AB007 | AB008 |
|---------------------------|-------|-------|
| Carbenicillin Sodium salt | 1g | 5g |

Storage conditions

Store at 2 - 8 °C

Stock solution: Prepare a stock solution with a concentration of 50 mg/ml in water and filter sterilize. Store the stock solution at -20°C. The working concentration varies from 20 - 100 µg/ml.

Application

- ✓ Used to select for ampicillin resistance in mutated and transformed cells (preferred over ampicillin due to its increased stability).
- ✓ Commonly used the elimination of Agrobacterium species after inoculation.

PRODUCT USE LIMITATION

This product is developed, designed and sold exclusively for research purposes and in vitro use only. The product was not tested for use in diagnostics or for drug development, nor is it suitable for administration to humans or animals. Please refer to www.canvaxbiotech.com for Material Safety Data Sheet of the product.



Tallaght Business Park Whitestown, Dublin 24, Ireland D24 RFK3

Tel: (01) 4523432 Fax: (01) 4523967 Web: www.labunlimited.com

Quatro House, Frimley Road, Camberley, United Kingdom GU16 7ER

Tel: 08452 30 40 30 Fax: 08452 30 50 30 E-mail: info@labunlimited.com E-mail: info@labunlimited.co.uk Web: www.labunlimited.co.uk





Notes



Tallaght Business Park Whitestown, Dublin 24, Ireland D24 RFK3

Tel: (01) 4523432 Web: www.labunlimited.com

Quatro House, Frimley Road, Camberley, United Kingdom GU16 7ER

Tel: 08452 30 40 30 Fax: (01) 4523967 Fax: 08452 30 50 30
E-mail: info@labunlimited.com E-mail: info@labunlimited.co.uk Web: www.labunlimited.co.uk

